

Table Water-curtain Spray Cabinet

BGD 228

Product Description

As the paint sprayed from the gun cannot completely cover on the work, so it produces great amount of poisonous coating mist when spraying the works. To keep the air not polluted and protect the workers' health, we produce a Table Water-curtain Spray Cabinet for small scale spraying in labs.

This spray cabinet applies the latest design plan, using the so-called 'negative pressure principle', the dental plate and arc plate produce strong air flow when working, and makes the water becomes eddy to wash the indrawn coating mist, the gas will be exhausted by the fan, and the paint residue left in the water. In addition, the whole spray cabinet is made of stainless steel and equipped with a high-pressure centrifugal fan, and it possesses a small footprint, is easy operate, safe, easy to clean. This spray cabinet is capable to directly splash the residual coating mist to the water pool or water-curtain, the processing efficiency up to over 90%. The smell and residual coating mist produced during the spraying will be filtered by the water curtain and exhausted outside the spraying room through the fan, so as to realize the cleansing of the spraying environment and the protection of people's health, as well as increasing the cleanliness of the works.



Technical Specification

- Coating mist collecting system: consists of stainless steel water-curtain plate, annular tank, water-curtain, and dash plate. Water-curtain plate, made of 1.5mm thickness stainless steel, facing towards operator. Water flows on its surface without break and pounding, maintaining a 2mm thickness water film. Most coating mist fully mixed with the water on the water curtain then flow into the annular tank, then filtered by the filter in the inlet of annual water pump.
- Water supply system: consists of annual water pump, valve, overflow channel and pipes.
- Exhausting system: Consists of Baffle-type steam separator, centrifugal exhaust fan, several exhaust pipe and fan holder, belonging to large flow and low thickness exhaust. Steam separator with maze structure fixed behind the water-curtain plate, capable to efficiently separate and condense the mist in the air, then flow back to the annual tank in case over lost fluid.

Table Water-curtain Spray Cabinet

BGD 228

Main Technical Parameters

- Material: Stainless Steel
- Overall Size: 810×750×1100 (L×W×H)
- Working Room Size: 600×500×380 (L×W×H)
- Exhaust Air Rate: 12m/s
- Fan: Single-phase centrifugal fan, power 370W
- Water Curtain Size: 600×400mm (L×W)
- Samples Holder Size: 595×200mm (L×W)
- Power Supply: 220V 50HZ
- Length of Air Duct: 2 m
- Ordering Information: BGD 228 - Table Water-curtain Spray Cabinet

Disclaimer

The information given in this sheet is not intended to be exhaustive and any person using the product for any purpose other than that specifically recommended in this sheet without first obtaining written confirmation from us as to the suitability of the product for the intended purpose does so at his own risk. Whilst we endeavour to ensure that all advice we give about the product (whether in this sheet or otherwise) is correct we have no control over either the quality or condition of the product or the many factors affecting the use and application of the product. Therefore, unless we specifically agree in writing to do so, we do not accept any liability whatsoever or howsoever arising for the performance of the product or for any loss or damage (other than death or personal injury resulting from our negligence) arising out of the use of the product. The information contained in this sheet is liable to modification from time to time in the light of experience and our policy of continuous product development