

Condensation Climatic Chambers

Cyclic Condensation-water Climatic Chamber BGD 876

Product Description

BGD 876 Cyclic Condensation-water Climatic Chamber is designed according to ISO 6270-2 (Paints and varnishes -- Determination of resistance to humidity -- Part 2: Condensation (in-cabinet exposure with heated water reservoir). This standard is different with Part 1"Continuous condensation", but defines three cyclic programs to simulate different humidity environments:

- CH: Condensation atmosphere with constant humidity
- AHT: Condensation atmosphere with alternating humidity and air temperature
- AT: Condensation atmosphere with alternating air temperature

It can give much better supplements and explanation for corrosion degree and defect types of coating used in different environments. In comparison to Part 1 - single-sided exposure this method adds a condensation atmosphere with alternating humidity and air temperature, and is mainly used to evaluate paints anti-corrosion ability under different environmental conditions.



Standards

ISO 6270-2 (paints and varnishes - Determination of resistance to humidity - Part 2: Condensation (in-cabinet exposure with heated water reservoir)



Condensation Climatic Chambers

Cyclic Condensation-water Climatic Chamber BGD 876

Technical Specification

Condensation Test Atmospheres

| Test Atmosphere | | | Cycle Duration | | Conditions in working chamber after reaching equilibrium | |
|---|--|------|---|-------|--|--|
| Туре | | Code | Test Period(s) | Total | Air Temperature | Relative Humidity |
| Constant Humidity condensation atmosphere | | СН | From warm-up to end of exposure | | 40±3°C | Approx 100% with condensation on test specimens |
| Alternating condensation atmosphere | With alternation of humidity and air temperature | AHT | 8h including warm-up | 24h | 40±3°C | Approx 100% with condensation on test specimens |
| | | | 16h including warm-up | | 18 to 28 °C | Approaching ambient |
| | With alternation of air temperature | AT | 8h including warm-up | 24h | 40±3°C | Approx. 100% with condensation on test specimens |
| | | | 16h including cooling down (chamber closed) | | 18 to 28°C | Approx. 100% (approximately saturated) |



Condensation Climatic Chambers

Cyclic Condensation-water Climatic Chamber BGD 876

Main Technical Parameters

- Designed with automatic ventilator, can control chamber ventilation automatically
- Operator can set three different working conditions: humidification, cooling with ventilation, cooling without ventilation and its corresponding temperature & running time. Any program including total test time and cycle times can be set freely
- Automatic water supply
- Protector for water temperature ceiling
- A glass observation window is designed in chamber, operator can observe specimens test status directly

Parameters:

- Test Specimens Capacity: 20 pcs (150mm×70mm standard size)
- Temperature Range of Working Room: RTD~60° (Dry or condensation)
- Overall Size: 800mm×750mm×1420mm (L×W×H)
- Weight: 180 kg
- Power: 220V 50HZ; Max. power of machine: 2.0KW
- Ordering Information: BGD 876 Cyclic Condensation-water Climatic Chamber

Disclaimer

The information given in this sheet is not intended to be exhaustive and any person using the product for any purpose other than that specifically recommended in this sheet without first obtaining written confirmation from us as to the suitability of the product for the intended purpose does so at his own risk. Whilst we endeavour to ensure that all advice we give about the product (whether in this sheet or otherwise) is correct we have no control over either the quality or condition of the product or the many factors affecting the use and application of the product. Therefore, unless we specifically agree in writing to do so, we do not accept any liability whatsoever or howsoever arising for the performance of the product or for any loss or damage (other than death or personal injury resulting from our negligence) arising out of the use of the product. The information contained in this sheet is liable to modification from time to time in the light of experience and our policy of continuous product development